


A scanning electron micrograph (SEM) showing a dense population of rod-shaped bacteria. The bacteria are oriented in various directions, some appearing as single cells and others as small clusters. Many of the rods have fine, hair-like flagella extending from their surfaces. The background is dark, making the light-colored bacterial structures stand out.

Counting Plates and Calculating Results

Counting colonies

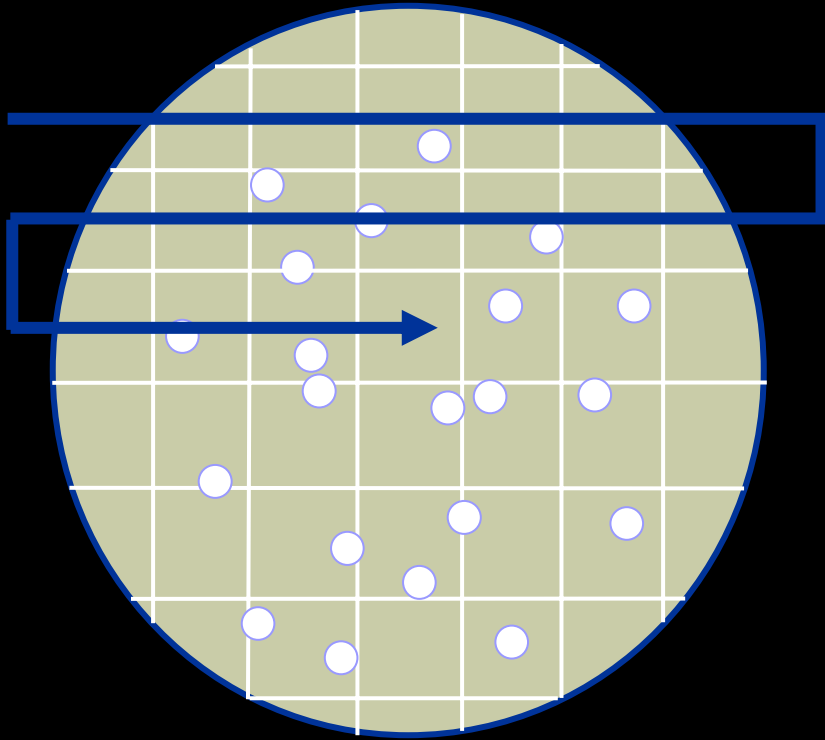

- Count colonies with a preset pattern using grid lines
- Recognize the difference between two or more colonies that have grown into contact with each other




Counting checks

- After one count, turn the plate 90 degrees and count again.
- Continue counting until counts agree within 5 percent.
- Have a second analyst check your count
- A count $>$ the ideal range: Estimate the count, if possible rather than “TNTC”



Ideal ranges of colony counts

Test type	Number of colonies
Fecal coliforms	20 – 60
Enterococci	20 – 60
Total coliforms	20 – 80
<i>Escherichia coli</i>	20 – 80
Fecal streptococci	20 – 100

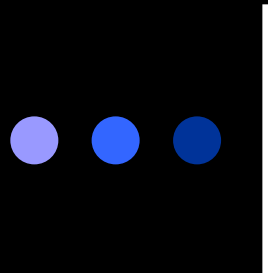

Basic formula

$$\text{Colonies / 100 mL} = \frac{\text{Number of colonies X 100}}{\text{Volume of sample filtered}}$$

Refer to Calculating Counts handout for different situations



Other scenarios

- Two or more filters within ideal range
 - Sum counts and sum volumes
Plate counts of 25 and 75 on 10 and 30 mL for *E. coli*:
 $(25+75)/(10+30)*100 = 250$ cfu/100 mL
- All filters above and/or below ideal range
 - Sum counts and volumes for all plates
Plate counts of 10 and 90 on 10 and 30 mL for *E. coli*:
 $(10+90)/(10+30)*100 = 250$ cfu/100 mL



Other scenarios

- No colonies

- Assume 1 colony on largest volume plated
- Code as “<” calculated concentration

No colonies on 3, 10, or 30 mL plates for *E. coli*:

$$(1)/(30)*100 = <3 \text{ cfu}/100 \text{ mL}$$

- All filters too numerous to count (TNTC)

- Assume maximum of ideal range on smallest dilution plated
- Code “>” calculated concentration

Lawn of growth on 3, 10, and 30 mL plates for *E. coli*:

$$(80)/(3)*100 = >2,700 \text{ cfu}/100 \text{ mL}$$

Reading the MPN table

# Large Wells Positive	IDEXX Quanti-Tray®/2000 MPN Table (per 100ml)																	
	# Small Wells Positive																	
	0	1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17
0	<1	1.0	2.0	3.0	4.0	5.0	6.0	7.0	8.0	9.0	10.0	11.0	12.0	13.0	14.1	15.1	16.1	17.1
1	1.0	2.0	3.0	4.0	5.0	6.0	7.1	8.1	9.1	10.1	11.1	12.1	13.2	14.2	15.2	16.2	17.3	18.3
2	2.0	3.0	4.1	5.1	6.1	7.1	8.1	9.2	10.2	11.2	12.2	13.3	14.3	15.4	16.4	17.4	18.5	19.5
3	3.1	4.1	5.1	6.1	7.2	8.2	9.2	10.3	11.3	12.4	13.4	14.5	15.5	16.5	17.6	18.6	19.7	20.8
4	4.1	5.2	6.2	7.2	8.3	9.3	10.4	11.4	12.5	13.5	14.6	15.6	16.7	17.8	18.8	19.9	21.0	22.0
5	5.2	6.3	7.3	8.4	9.4	10.5	11.5	12.6	13.7	14.7	15.8	16.9	17.9	19.0	20.1	21.2	22.2	23.3
6	6.3	7.4	8.4	9.5	10.6	11.6	12.7	13.8	14.9	16.0	17.0	18.1	19.2	20.3	21.4	22.5	23.6	24.7
7	7.5	8.5	9.6	10.7	11.8	12.8	13.9	15.0	16.1	17.2	18.3	19.4	20.5	21.6	22.7	23.8	24.9	26.0
8	8.6	9.7	10.8	11.9	13.0	14.1	15.2	16.3	17.4	18.5	19.6	20.7	21.8	22.9	24.1	25.2	26.3	27.4
9	9.8	10.9	12.0	13.1	14.2	15.3	16.4	17.6	18.7	19.8	20.9	22.0	23.2	24.3	25.4	26.6	27.7	28.9
10	11.0	12.1	13.2	14.4	15.5	16.6	17.7	18.9	20.0	21.1	22.3	23.4	24.6	25.7	26.9	28.0	29.2	30.3
11	12.2	13.4	14.5	15.6	16.8	17.9	19.1	20.2	21.4	22.5	23.7	24.8	26.0	27.2	28.3	29.5	30.7	31.9
12	13.5	14.6	15.8	16.9	18.1	19.3	20.4	21.6	22.8	23.9	25.1	26.3	27.5	28.6	29.8	31.0	32.2	33.4
13	14.8	16.0	17.1	18.3	19.5	20.6	21.8	23.0	24.2	25.4	26.6	27.8	29.0	30.2	31.4	32.6	33.8	35.0
14	16.1	17.3	18.5	19.7	20.9	22.1	23.3	24.5	25.7	26.9	28.1	29.3	30.5	31.7	33.0	34.2	35.4	36.7
15	17.5	18.7	19.9	21.1	22.3	23.5	24.7	25.9	27.2	28.4	29.6	30.9	32.1	33.3	34.6	35.8	37.1	38.4
16	18.9	20.1	21.3	22.6	23.8	25.0	26.2	27.5	28.7	30.0	31.2	32.5	33.7	35.0	36.3	37.5	38.8	40.1
17	20.3	21.6	22.8	24.1	25.3	26.6	27.8	29.1	30.3	31.6	32.9	34.1	35.4	36.7	38.0	39.3	40.6	41.9
18	21.8	23.1	24.3	25.6	26.9	28.1	29.4	30.7	32.0	33.3	34.6	35.9	37.2	38.5	39.8	41.1	42.4	43.8
19	23.3	24.6	25.9	27.2	28.5	29.8	31.1	32.4	33.7	35.0	36.3	37.6	39.0	40.3	41.6	43.0	44.3	45.7
20	24.9	26.2	27.5	28.8	30.1	31.5	32.8	34.1	35.4	36.8	38.1	39.5	40.8	42.2	43.6	44.9	46.3	47.7
21	26.5	27.9	29.2	30.5	31.8	33.2	34.5	35.9	37.3	38.6	40.0	41.4	42.8	44.1	45.5	46.9	48.4	49.8
22	28.2	29.5	30.9	32.3	33.6	35.0	36.4	37.7	39.1	40.5	41.9	43.3	44.8	46.2	47.6	49.0	50.5	51.9
23	29.9	31.3	32.7	34.1	35.5	36.8	38.3	39.7	41.1	42.5	43.9	45.4	46.8	48.3	49.7	51.2	52.7	54.2
24	31.7	33.1	34.5	35.9	37.3	38.8	40.1	41.7	43.1	44.6	46.0	47.5	49.0	50.5	52.0	53.5	55.0	56.5
25	33.6	35.0	36.4	37.9	39.3	40.8	42.2	43.7	45.2	46.7	48.2	49.7	51.2	52.7	54.3	55.8	57.3	58.9
26	35.5	36.9	38.4	39.9	41.4	42.8	44.3	45.9	47.4	48.9	50.4	52.0	53.5	55.1	56.7	58.2	59.8	61.4
27	37.4	38.9	40.4	42.0	43.5	45.0	46.5	48.1	49.6	51.2	52.8	54.4	56.0	57.6	59.2	60.8	62.4	64.1
28	39.5	41.0	42.6	44.1	45.7	47.3	48.8	50.4	52.0	53.6	55.2	56.9	58.5	60.2	61.8	63.5	65.2	66.9
29	41.7	43.2	44.8	46.4	48.0	49.6	51.2	52.8	54.5	56.1	57.8	59.5	61.2	62.9	64.6	66.3	68.0	69.8
30	43.9	45.5	47.1	48.7	50.4	52.0	53.7	55.4	57.1	58.8	60.5	62.2	64.0	65.7	67.5	69.2	71.0	72.8